LUMATEC

---

# SUPERLITE M 03

---

FORENSIC  
MULTISPECTRAL  
LED STRIP LIGHT

Operating Instructions

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[EN] These operating instructions belong with this product. They contain important information for putting it into service and operating it. This should be noted also when this product is passed on to a third party.

Therefore look after these operating instructions for future reference!

A list of contents with the corresponding paged numbers can be found in the index on page 1.

LUMATEC prepared these instructions for experienced forensic scientists, process engineers, technicians and manufacturing personnel. If you are new to light applications, do not understand the instructions or have any further questions and concerns, you can find our contact information on the last page of this booklet.

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#### Part No

2500.0000	SUPERLITE M 03	[LED strip light]
2500.0010	SUPERLITE M 03	[LED strip light with case]
9200.0139	SUPERLITE M 03	[Telescopic support rod with ball head / spare part]
9200.0140	SUPERLITE M 03	[Charge for LED strip light / spare part]

---

**TABLE OF CONTENTS**

1. Intended Use..... 2

2. Safety Instructions..... 2

3. Equipment Overview ..... 2

4. Equipment Installation ..... 4

5. Operation ..... 4

6. Accessories and Spare Parts ..... 6

7. Technical Data..... 7

---

## 1. INTENDED USE

The SUPERLITE **M** 03 LED Strip Light is solely intended for visualizing of shape traces on surfaces.

---

## 2. SAFETY INSTRUCTIONS

Before you start working with the SUPERLITE **M** 03, please read these operating instructions completely.

### Power Supply

For charging and power supply in mains operation, only use the power supply unit / charger included in the delivery. This automatically adjusts to all AC currents from 110 to 240 Volts and 50 to 60 Hz.

### Repairs

Do not attempt to modify or repair the product. All repairs, including replacing the permanently installed rechargeable battery, may only be carried out by authorized personnel.

### Explosive Environments

This device is not meant for operation in explosive environments.

### Humid Environments

The strip light is dust-tight and protected against splash water from all directions (IP 64), but must not be immersed in water.

### Protection against Strong Optical Radiation

Do not look directly into the strip light. Do not direct the light beam into the eyes of humans or animals. If someone is still blinded by the glare, medical help should be requested immediately.

## Danger of Fire

Do not cover or pack the device during operation to avoid excessive heat build-up. Switch the device off during transport.

---

## 3. EQUIPMENT OVERVIEW

When strip light falls flat on a surface, even the smallest structures become visible through high-contrast shadows. The brightness, color and color saturation of the grazing light can be individually adjusted and thus optimally adapted to the surface. There are 5 memory spaces available for saving your own settings. With the help of the telescopic support rod (length 21 to 91 cm) combined with an extremely strong ball head, the light can be easily guided and placed on the floor or a wall. Tiny feet prevent traces from being smeared even when the unit is put down on the ground.

Alternatively, there are 1/4" UNC threaded holes in the middle on the top and bottom for mounting on a tripod.

The light is equipped with an integrated Li-Ion battery for mobile use of up to 120 minutes at 50 % brightness. The scope of delivery includes the required charger with USB-C socket, which is also used for stationary mains operation.

SUPERLITE M 03 LED Strip Light

Application: Print Shoe Size: EU 43 / UK 9

Control Panel:

Display

Button Color

Button Saturation

Button Brightness

Button User Setting

Control Panel

USB-C Charging Socket

On/Off Rocker Switch

Telescopic Support Rod

Tiny Feet (8x)

Ball Head

Clamping Lever

---

## 4. EQUIPMENT INSTALLATION

### Assembly of the Telescopic Support Rod

Open the clamp on the ball head using the lateral clamp lever so that the ball can move freely with the connected 1/4" tripod screw with plate. Turn the screw clockwise into the silver threaded hole on top of the housing until it stops and bring the support rod into a vertical position. Close the clamp lever on the ball head again.

### Adjust the Length of the Tele- scopic Support Rod

The length of the support rod can be extended or shortened by pulling and pushing. When the rod is in a vertical position, hold the ball head with one hand while you adjust the length at the other end with the other hand.

### Angle Adjustment

To adjust the angle, open the clamp on the ball head using the lateral clamp lever, set any angle and close the clamp lever again.

Please note that if the rod is fully extended and not in a vertical position, the setup may tip over when put down later.

For an extremely angled position, ideally with a short length, move the rod in the direction of the side recess of the ball head when adjusting the angle.

### Mounting on a Tripod or Tripod Head

Instead of the telescopic support rod, the strip light can also be mounted on

a tripod or tripod head using the silver 1/4" threaded hole on the top and bottom.

Note: Do not mount the strip light at eye level so that people are not blinded, but rather close to the surface that is to be examined (e.g. floor or wall)!

### Dismantling the Support Rod or Tripod

Dismantling is carried out in the reverse order of assembly, see above.

---

## 5. OPERATION

### Switching On and Off

The rocker switch, which is located at the right outside of the housing is used to turn the strip light on and off, position 1 and 0. A protective cover on the side prevents it from being switched on accidentally, e.g. during transport. After switching on, the current color, brightness (with dim: 25%) and battery level (battery symbol) are shown on the display.

### Charging

Next to the rocker switch is the USB-C charging socket.

Switch the device off (rocker switch in position 0) and plug the USB-C cable into the unit's charging socket on one side and into the USB-C charger on the other side. Then plug the USB-C charger into a power outlet. During charging, the current battery level is shown in % on the display. The charging process from 5-100 % takes about 100-120 minutes. When the charging process is complete, "Charging Done" appears on the display.

Then disconnect the USB-C cable from the unit.

Note: Only use the original charger included in delivery!

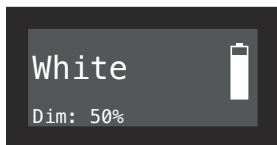
## Mains Operation

Connect the device to the USB-C charger as described under “Charging” and switch the device on using the lateral rocker switch (position 1). The display then switches to operating mode.

## Battery Operation

After charging, disconnect the USB-C cable from the unit and switch the device on (rocker switch in position 1). The display then switches to operating mode.

## Display

The current color, saturation or brightness value is

shown on the display depending on the last button press. In addition, the current brightness value is always shown in % at the bottom left. The battery status display is shown on the right edge.

## Operating Concept of the Color, Color Saturation and Brightness Buttons

By briefly pressing the respective button, you change the settings in large steps, e.g. for color saturation and brightness in 25 % steps.

By pressing the button for a longer time, that means pressing and holding it, you change the value in small steps, for example 1 %.

By pressing and holding it again, you change the direction of the setting, e.g. the values displayed will no longer be greater but smaller.

## The Color Button

By pressing and holding, color values between 0 and 360 can be set.

By briefly pressing, you can switch between the current color value (shown in the display with “Color:” 0-360) and white (shown in the display with “White”). This allows you to quickly assess whether the current color value (0-360) leads to a better result than white.

Tip: When setting the color, it is advisable to set the color saturation (button to the right of color) to 100 % so that changes can be clearly seen.

## The Saturation Button

A short press on this button reduces the color saturation by 25 % in one step.

A long press changes the color saturation in 1 % increments. An additional long press changes the direction of the setting (increasing or decreasing values)

The current % value is shown in the display after “Sat:”.

Note: If the color is set to “White”, the color saturation cannot be adjusted because the color saturation of white is 0 by definition. Press the color button to select a “colorful” color and then press the color saturation button again.

## The Brightness Button

A short press on this button reduces the brightness by 25 % in one step.

A long press changes the brightness in 1 % increments. Pressing and holding again changes the direction of the setting (increasing or decreasing values)

The current % value is shown in the display after “Dim:”.

## The User Setting Button

By pressing briefly this button, the 5 previously saved, user settings are called up one after the other. “CUS:” and the corresponding number appear in the display.

To save a new user setting, select the memory space you want to overwrite, for example CUS: 5.

Tip: The memory space you are overwriting is now shown in the bottom right of the display with “CUS:” and the corresponding number.

Now make the desired settings using the color, brightness and color saturation buttons. Then press and hold the user setting button until ...SAVED... appears in the display.

## 6. ACCESSORIES AND SPARE PARTSE

9200.0138	Case for SUPERLITE <b>M</b> 03
9200.0139	Telescopic support rod / length 25 to 95 cm / spare part
9200.0140	Charger for SUPERLITE <b>M</b> 03 / USB-C socket / spare part

Image above:

Extended SUPERLITE **M** 03 set with  
case 9200.0138 and other devices from  
the SUPERLITE **M** 05 system

---

## TECHNICAL DATA

Model:	SUPERLITE <b>M</b> 03 / Part No. 2500.0000
Luminous flux::	max. 1656 lm (measured at color setting "White")
Optical power:	max. 9,1 W (measured at color setting "White")
Dimensions: (L / W / H):	410 mm / 67 mm / 28 mm
Weight:	670 g
Battery:	14.4 V, 3.5 Ah, Lithium-Ion
Number of selectable colors:	360
Adjustable saturation settings:	0 % to 100 % (in 1 % or 25 % steps)
Adjustable brightness settings:	0 % to 100 % (in 1 % or 25 % steps)
Custom defined memory spaces:	5
Battery runtime:	120 min (at 50 % brightness)
Operating temperature:	+ 10 °C to + 35 °C
Storage temperature:	- 10 °C to + 60 °C
Relative humidity:	< 75 % (< 30 °C) < 50 % (> 30 °C)
Input power supply / charger:	100-240 VAC, 50 - 60 Hz
Output power supply / charger:	max. 65 W PD / USB-C
Conformity:	EN 55015, EN 60598-1, EN 62471 Low Voltage Directive 2014/35/EU Electromagnetic Compatibility 2014/30/EU Ecodesign Directive 2009/125/EC RoHS Directive 2011/65/EU

### Declaration of conformity

This unit conforms to all applicable EC Directives and corresponding harmonised standards. A written declaration of conformity can be supplied on request.

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LUMATEC GmbH